


- Remove plastic wrapping.
- Before opening, tap the base of the box on the bench to ensure all tips are fully nested.

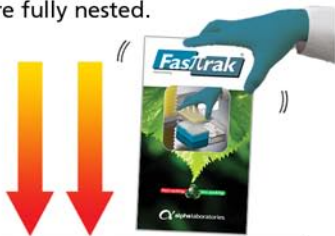

1.

- Undo the paper tab and open the box by lowering both sides simultaneously.
- Grasp transfer top. Lift tips straight up and clear of stack.



2.

- Align transfer top over rack and place onto rack.



3.

- Push down on transfer top to release refill plate and click onto rack. Return transfer top to Fastrak[®] refill stack.




4.

- Close the Fastrak[®] carton, lifting both sides simultaneously and re-seal the paper tab.
- Once in use, do not invert the pack.



5.

- Store boxes upright to ensure good tip nesting.

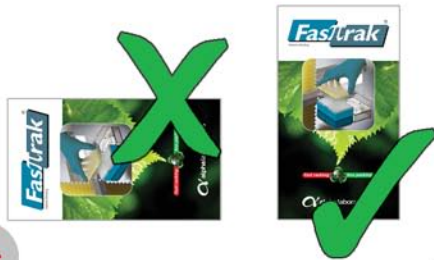

6.